

How to print the column as you want in calorimeter software

1. Click "View History" to check the results you analyzed as Figure 1;

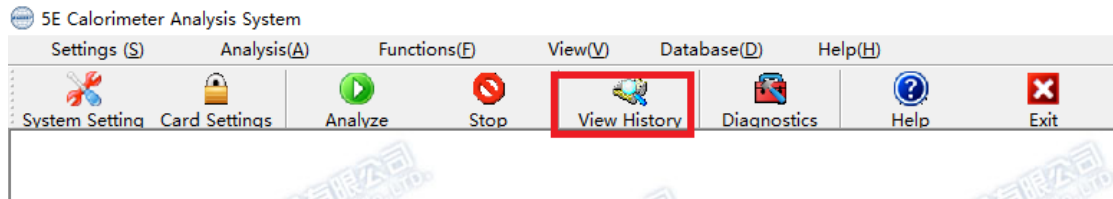


Figure 1

2. Right click you mouse, you will get choice as showed in Figure 2, choose "Datagrid Setting";

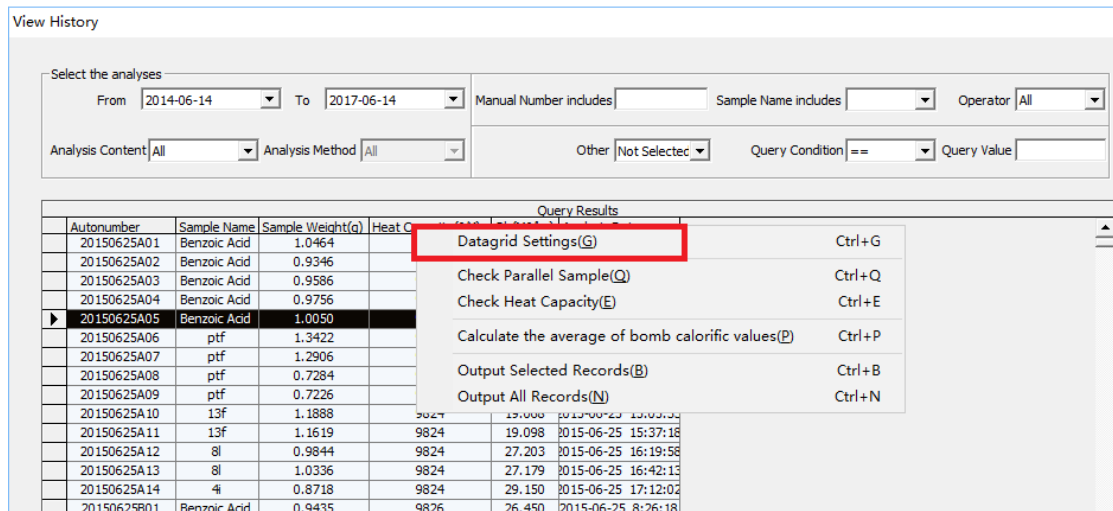


Figure 2

3. Then choose the column you want to display or hide as in Figure 3, and click ok to save;

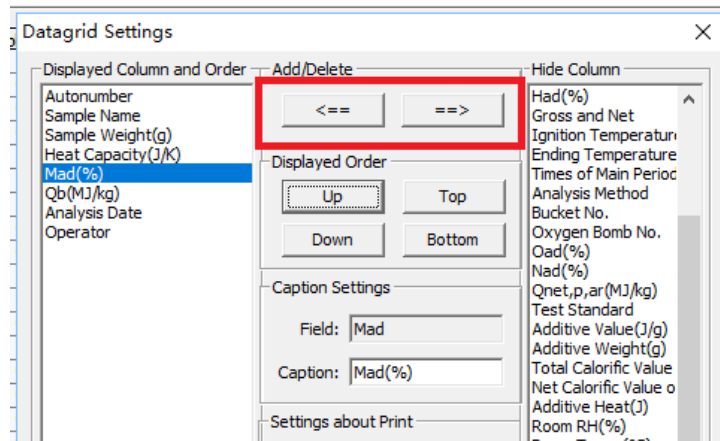


Figure 3

4. Select the records that you want to print, then click "Print the Selected Records" as in Figure 4

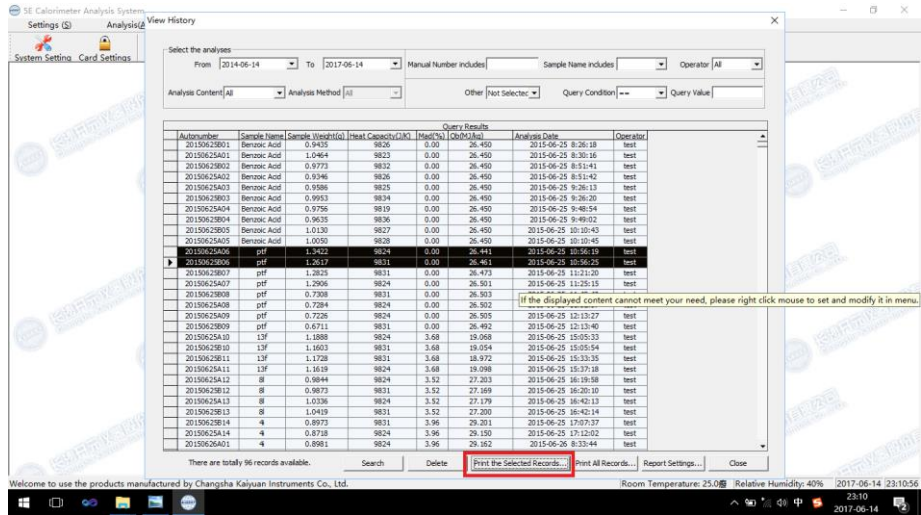


Figure 4

4. The columns will be printed as you set as showed in Figure 5

Analysis Data Report

Equipment Model: 5E-C5500				Equipment Serial:			
Autonumber	Sample Name	Sample Weight (g)	Heat Capacity (J/K)	Mad (%)	Qb (MJ/kg)	Analysis Date	Operator
20150625A06	ptf	1.3422	9824	0.00	26.441	2015-06-25	test
20150625B06	ptf	1.2617	9831	0.00	26.461	2015-06-25	test

Company:

Operator:

Auditor:

Figure 5